


ReliChrom™

Ready to use
pre-packed columns

ReliChrom™

Ready to use pre-packed columns

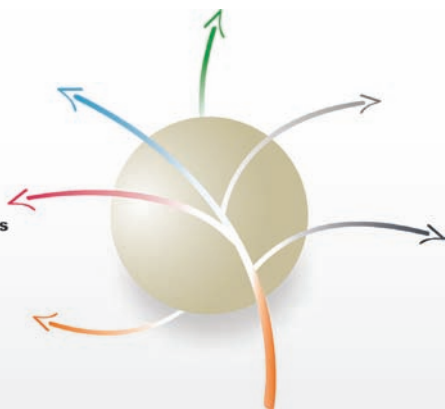

Resindion S.r.l.

A Subsidiary of  MITSUBISHI CHEMICAL

www.resindion.com

ReliChrom™

Ready to use pre-packed columns

Typical features


ReliChrom™ pre-packed columns have been developed by Resindion to provide a practical and reliable tool to screen **ReliSorb™ SS** highly porous polymeric matrices in chromatographic bioseparations.

ReliChrom™ columns have been conceived for a possible direct connection to almost all standard LC and HPLC systems through UNF 10-32 male (1/16" male) fittings. Each column contains 5 ml of packed resin and the resin bed height is 100 mm. These dimensions represent an excellent solution for the development of optimized purification methods.

ReliChrom™ column hardware is designed for a max operating pressure of 30 bar (430 psi) and for a 5 - 60 °C temperature range in operation. Overall column dimensions are 11.5 mm (8 mm ID) x 135 mm (100 mm HR); each column is equipped with 17 µm frits.

All the components of **ReliChrom™** columns are made of polymeric materials: polypropylene (PP) and high density polyethylene (HD-PE).

ReliChrom™ columns are chemically stable in all pH range (1 - 14), in high salt concentration buffer and in common solvents (avoid use of strong oxidizing agent and halogenated solvents). Proper plug-stoppers are used to grant the packed media stability during storage time.

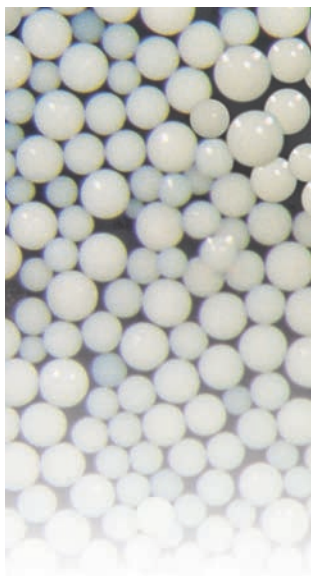

SALES
ON
THE
Resindion
Shop

ReliChrom™ columns characteristics

ReliSorb™ SS* packed volume:	5.0 ml
Resin bed depth:	100 mm
ReliSorb™ SS particle size range:	50 - 150 µm
Operating pH range:	1 - 14
Operating temperature range:	5 - 60 °C
Storage temperature range:	5 - 30 °C (4 - 6 °C once used)

* **ReliSorb™** are also available in bulk in the following particle size: SS (50-150 µm), standard (75-200 µm) and EB (200-500 µm).



Characteristics of ReliChrom™ packed columns

NAME	FUNCTIONAL GROUP	IONIC FORM AT DELIVERY	Min. ION EXCHANGE CAPACITY (meq/ml)	BASIC APPLICATION	DYNAMIC BINDING CAPACITY (DBC)		
					BSA (mg/ml)	Lys ⁵ (mg/ml)	Papaine ^d (mg/ml)
CM400/SS	Carboxyl	H ⁺	0.15	CEX	—	min. 30	—
SP400/SS	Sulphopropyl	Na ⁺	0.10	CEX	—	min. 40	—
DA400/SS	Tertiary amine	free base	0.30	AEX	min. 30 ^a	—	—
QA400/SS	Quaternary ammonium	Cl ⁻	0.30	AEX	n. a.	—	—
IDA400/SS	Iminodiacetic	Ni ²⁺	100 (μmol/ml Ni ²⁺) ¹	IMAC	—	—	10
BU400/SS	Butyl	—	—	HIC	min. 30 ^b	—	—
PH400/SS	Phenyl	—	—	HIC	min. 30 ^b	—	—
OD400/SS	Octadecyl	—	—	HIC	min. 20 ^b	—	—

¹Ni²⁺ ionic form is applied only for ReliChrom™ pre-packed columns; different ionic forms are also available upon request.

^aFeed solution: 10 g/l BSA in 20 mM TRIS - HCl buffer, pH 7; flow rate = 150 cm/h

^bFeed solution: 10 g/l BSA in 20 mM phosphate buffer, pH 7 + (NH₄)₂SO₄ 2 M; flow rate = 150 cm/h

^cFeed solution: 8 g/l Lysozyme in 20 mM sodium acetate buffer, pH 5; flow rate = 150 cm/h


^dFeed solution: 20 g/l Papaine crude extract in 20 mM phosphate buffer, pH 7.2 + NaCl 200 mM; flow rate = 150 cm/h

Supply conditions:

- IEX and IMAC columns: EtOH 20% aqueous solution + NaCl (final concentration 150 mM)

- HIC columns: EtOH 20% aqueous solution

Pressure drop in water at 25 °C

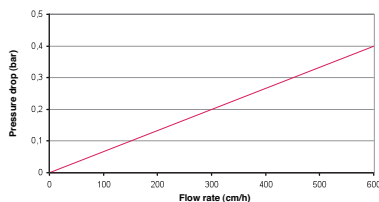

Column dimensions:

i.d. = 0.8 cm
length = 10 cm
area = 0.5 cm²
ReliSorb™ SS packed volume = 5 ml

Handling and storage

- Solutions to be treated, eluents, regenerant and storage solutions should always be freshly prepared, filtered (0.45 μm membrane filter) and degassed.
- Prior to connecting the column to any liquid supply, remove air bubbles from the inlet tube.
- Before the column first use it is recommended to measure, at various flow rates, the pressure drop on the ReliChrom™ induced by the chosen eluent, in order to obtain reference data.
- Set the maximum linear flow rate to a value appropriate to the resin considered.
- A significant pressure drop increase after repeated cycles and under the same conditions could indicate the presence of impurities: in this case the column has to be submitted to standard regeneration or cleaning in place procedures.
- After the operation and in case of long storage, the regeneration step has to be followed by a resin conditioning with storage solution. After use, ReliChrom™ columns have to be stored at 4 – 6 °C.

ReliChrom™

Visit our
web site
and
discover
the
Resindion
Shop

Material Support File



Resindion S.R.L. is an UNI EN ISO 9001:2008 certified Company. Technical and basic regulatory documents, supplied with each **ReliChrom™** delivery, are related to the Handling Instructions and Column Performance Report (HICPR).


Resindion S.R.L. can supply the information related to the **ReliSorb™ SS** packed resins (i.e. Technical Data Sheet, Certificate of Analysis, Material Safety Data Sheet).

Customer service

Resindion S.R.L. qualified technical and marketing team grants all the necessary customer assistance on **ReliChrom™** selection and application.

Please consult Resindion at: customerservice@resindion.com or technicalservice@resindion.com

Resindion Product Lines

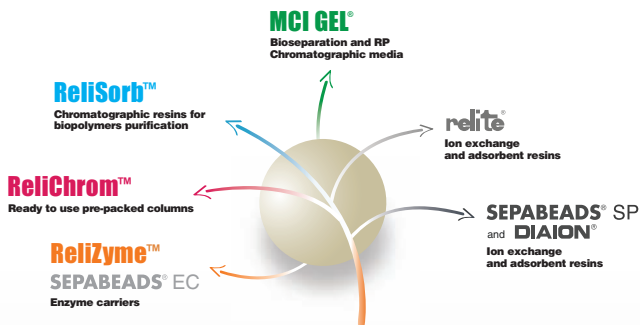

Resindion S.r.l.

A Subsidiary of MITSUBISHI CHEMICAL

Resindion S.R.L.

Via Roma 55
20082 Binasco (MI), Italy
Phone: +39 02 9001301
Fax: +39 02 90090774
E-mail: info@resindion.com
Web site: www.resindion.com

Mitsubishi Chemical Corporation

Palace Building
1-1, Marunouchi 1-chome, Chiyoda-ku
Tokyo 100-8251, Japan
Phone: +81 3 6748 7146
Fax: +81 3 3286 1209
Web site: www.diaion.com



Neither this publication, nor any product in this brochure shall be construed as recommending the infringement of any patent, nor extending any license, expressed or implied, nor assuming any liability under any issued or pending patent. The data presented have been carefully compiled from Resindion records which are believed to be accurate and reliable. However, no warranties or representations with respect hereto are made, nor is freedom from any patent to be inferred.